



10451 ROSELLE STREET, #300, SAN DIEGO, CA 92121
TELEPHONE (858) 642-1988 • FAX (858) 642-1989
WWW.ATSBIO.COM • ATS@ATSBIO.COM

Goat IgG-SAP
NON-TARGETED SAPORIN CONTROL MOLECULE

[pre-immune goat IgG antibody]-saporin

Catalog Number: IT-19
Quantity: 25 micrograms, 100 micrograms, 250 micrograms
Format: PBS (0.14 M Sodium Chloride; 0.003 M Potassium Chloride; 0.002 M Potassium Phosphate; 0.01 M Sodium Phosphate; pH 7.4), no preservative. Sterile-filtered.
Host: Goat

Background:

Controls are a vital part of the scientific procedure; without them it is difficult to isolate the specific effects from the non-specific or artifactual. This control molecule is the same molecular weight, consists of similar, comparable materials and is synthesized with the same protocols as the targeted immunotoxins. The difference is the cell-specific targeting agents are replaced with "blanks," antibodies or peptides that have no specificity, and no ability to target cells. In short, they are the perfect control molecules for behavioral experiments with Advanced Targeting Systems' immunotoxins.

Specificity and Preparation:

This control conjugate (molecular weight 210 kDa) has no known specificity; it may react with cells that possess Fc receptors. Goat IgG-SAP is a chemical conjugate of pre-immune goat IgG antibody and the ribosome inactivating protein, saporin. The product is routinely tested by cytotoxicity assay.

Usage and Storage:


Goat IgG-SAP serves as a control for goat IgG-containing immunolesioning agents (Mab-ZAP, Rab-ZAP, Hum-ZAP, Rat-ZAP, Anti-M-ZAP). There may be lot-to-lot variation in material; working dilutions must be determined by end user.

Centrifuge material at low speed in microfuge to ensure all of solution is at bottom of tube. Vortex gently. The material should be stored at -20°C in undiluted aliquots. Material should be aliquoted to a convenient volume and quantity to avoid repeated freezing and thawing that can damage the protein content. Under these conditions, the material has a very stable shelf-life. Thawing should be done at room temperature or on ice. The thawed solution should remain on ice until use. The material can be handled safely using normal laboratory precautions.

Do not use a reducing agent (such as dithiothreitol, beta-mercaptoethanol or ascorbic acid) with this material. It will inactivate the toxin.

For disposal: autoclave, or expose to 0.2 M NaOH, materials that come into contact with the toxin.

Goat IgG-SAP
NON-TARGETED SAPORIN CONTROL MOLECULE

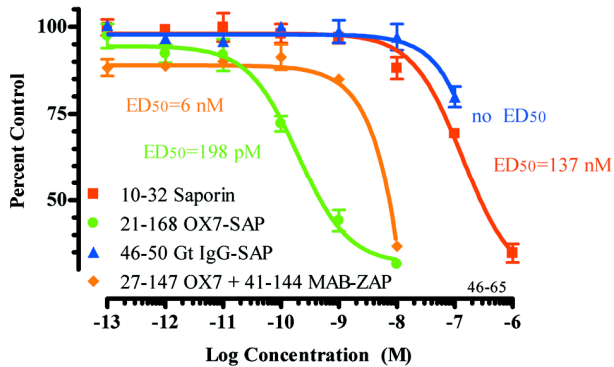

PC12 cells, a rat tumor cell line that expresses the Thy 1 receptor, were plated at 5000 cells per well and allowed to acclimatize overnight. The cells were exposed to OX7-SAP, OX7 Ab (diluted in 50 ng/well MAB-ZAP), and Goat IgG-SAP at the indicated concentrations for 72 hours. MTS/PMS (Promega) mixture was added and plates were read on a Molecular Diagnostics Spectramax plate reader with SoftMax software. Data was analyzed by Prism 4.0 software.

References:

1. Tazzari PL, Bolognesi A, De Toter D, Falini B, Lemoli RM, Soria MR, Pileri S, Gobbi M, Stein H, Flenghi L, *et al.* (1992) Ber-H2 (anti-CD30)-saporin immunotoxin: a new tool for the treatment of Hodgkin's disease and CD30+ lymphoma: *in vitro* evaluation. *Brit J Haemat* 81:203-211.
2. Dinota A, Tazzari PL, Michieli M, Visani G, Gobbi M, Bontadini A, Tassi C, Fanin R, Damiani D, Grandi M, *et al.* (1990) *In vitro* bone marrow purging of multidrug-resistant cells with a mouse monoclonal antibody directed against Mr 170,000 glycoprotein and a saporin-conjugated anti-mouse antibody. *Cancer Res* 50:4291-4294.
3. Thorpe PE, Brown AN, Bremner JA Jr, Foxwell BM, Stirpe F (1985) An immunotoxin composed of monoclonal anti-Thy 1.1 antibody and a ribosome-inactivating protein from *Saponaria officinalis*: potent antitumor effects *in vitro* and *in vivo*. *J Natl Cancer Inst* 75(1):151-159.

Safety:

Good laboratory technique must be employed for safe handling of this product.

This requires observation of the following practices:

1. Wear appropriate laboratory attire, including lab coat, gloves and safety glasses.
2. Do not pipet by mouth, inhale, ingest or allow product to come into contact with open wounds. Wash thoroughly any part of the body which comes into contact with the product.
3. Avoid accidental autoinjection by exercising extreme care when handling in conjunction with any injection device.
4. This product is intended for research use by qualified personnel only. It is not intended for use in humans or as a diagnostic agent. Advanced Targeting Systems is not liable for any damages resulting from the misuse or handling of this product.